VWR® ELECTRONIC PIPETTE

QUICK START GUIDE

EN / FR / DE / ES / IT / PT

For more information, please consult the full manual

VWR® ELECTRONIC PIPETTE

ENGLISH

Box contents

- VWR® Electronic pipette,
- Power Supply 5V with AC adapter and cable,
- 4 Battery side tags,
- Lubricant (except for 10 μ L and the multichannel models),
- Tip ejector extension (10 μ L model only),
- Quick Start Guide,
- Certificate of conformity (including bar-code sticker),
- 2-year warranty.

Installation

Prior to use, fully charge the VWR Electronic pipette; this will take approximately 3 hours. Charge by using one of the following:

- Power supply, included in box
- Stand adapter, for use with VWR Carrousel Pipette Stand, sold separately.

Note: Battery wake-up can take a few minutes.


Only use the original power transformer supplied by VWR. Use of an incompatible power supply can damage your VWR Electronic pipette.

VWR Electronic pipette uses a Lithium-Ion battery. Operation on the battery is done at user's risk only. Opening the pipette voids the warranty.

Getting started with VWR Electronic Pipette

VWR Electronic pipette has only two buttons to enable the pipette and reach all menu options:

- **Push-button:** set the volume by turning the button left or right and pipette with a simple push.
- **Mode-button:** access the pipette modes and readjustment functions.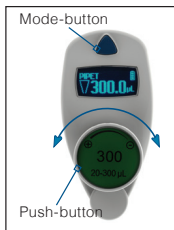

Using the push-button


To wake, or turn on the pipette, press the push-button.

The start-up screen appears.


The pipette will do a self-calibration test, the firmware resets the piston.

VWR Electronic pipette is preset on PIPET mode also known as standard pipetting.

Volume range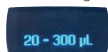

Version of firmware

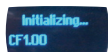Note: VWR Electronic pipette goes into sleep mode after 3 minutes of inactivity. The display will turn off. Simply press the push-button to reactivate the pipette. Your last setting and pipetting mode will appear on the screen.

To adjust the volume setting

- Quickly turn the push-button halfway, the screen will display "Click to change volume".
- Press the push-button; the volume on display will start blinking.
- You can now adjust the volume by turning the push-button either clockwise to decrease volume or counter clockwise to increase volume.
- Press the push-button once when finished to lock the volume.

To aspirate and dispense

- After the volume setting, just press the push-button to aspirate and press again to dispense.
- If you keep your thumb pressing the push-button until the end of the dispense cycle: the piston reset takes place after releasing the push-button again. It enables you to remove the tip from the liquid, without aspirating anything.

Using the mode-button

You can access the menu (pipetting modes, speed, configuration, quit) by pushing the mode-button. To choose one of the modes, turn the push-button and confirm your selection with a simple click!

The speed is set by default on speed 6 (maximum speed) but you can change aspiration and dispensing speed independently by reaching the speed menu.

Available modes:

- **PIPET:** standard pipette mode; the pipette aspirates and dispenses the volume set.
- **REPETITIVE:** pipette distributes a same volume repeatedly in predefined number of equal aliquots.
- **MIX:** standard pipette mode followed by a mixing phase (=repeatedly aspirating and dispensing), and optional standard pipette phase.
- **REVERSE:** a preset amount of liquid is added during aspiration. After delivery, the excess volume remains in the tip and should be discarded.

VWR® ELECTRONIC PIPETTE

- **CUSTOM:** allows to build a customized protocol to be transferred to the pipette using VWR protocol builder software. To download, go to vwr.com/e-pipette

Switching off the pipette

To switch off VWR Electronic pipette, press the push-button and hold for at least 5 seconds or it will switch off after 3 minutes of inactivity.

Other references

The Quick Start Guide provides you with elementary information to get started with the VWR Electronic pipette. Make sure you have read the Users' Guide, downloadable at: vwr.com/e-pipette for detailed information.

Visit vwr.com for additional information.

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LT801558/D © 12/2021
United States: VWR International LLC - 100 Matsonford Rd - Radnor, PA 19087 -