


Short user's manual

- VWR[®] UHP line
- VWR[®] EHP line
- VWR[®] Standard line
- VWR[®] EHP, fixed volume



This short user's manual is intended to provide a brief overview of the main features and operation of all VWR[®] Single- and Multi-channel Pipettors.

For detailed information, please refer to the User Manuals that can be found at www.vwr.com in several languages.

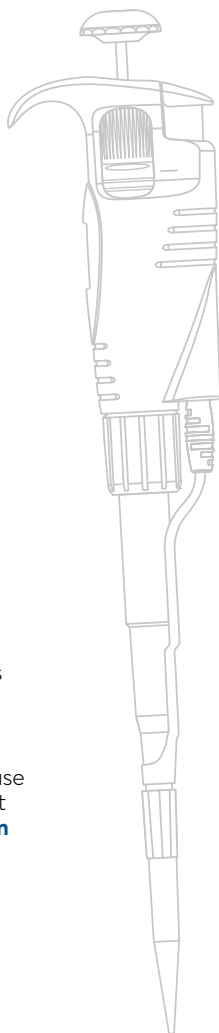


Getting started

Intended Use


VWR® Pipettors are volumetric instruments designed to measure and transfer liquids precisely and safely with the use of disposable pipet tips. The pipettors operate on the air-cushion principle (i.e., the aspirated liquid does not come in contact with the shaft or plunger of the pipettor).

They can accommodate volumes from 0.1 µL to 10,000 µL depending on the model. They are available in single-, 8- and 12-channel variable volume as well as fixed volume single-channel versions.

Variable volume pipettors

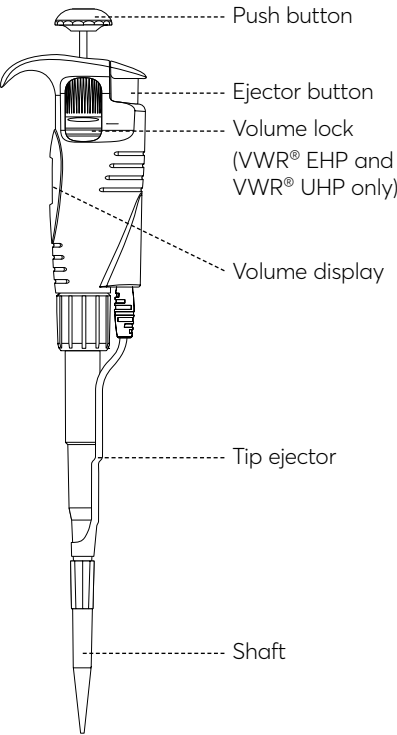


VWR® EHP, fixed volume

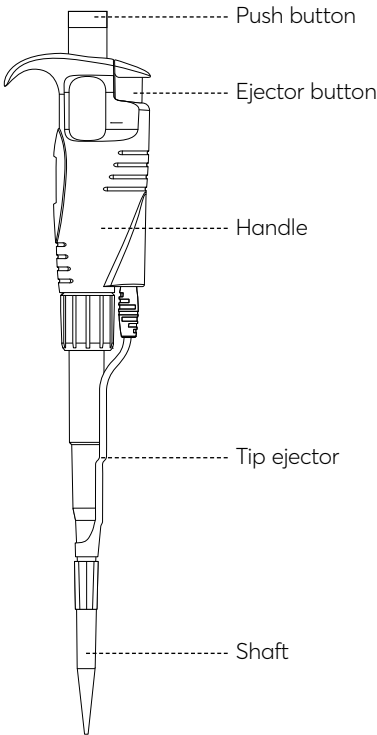

Step 1

Set the volume

■ VWR® EHP ■ VWR® Standard

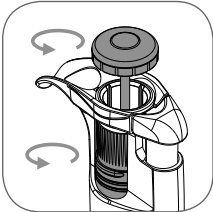

2	1
1	2
6	5

2.16 μ L 1.25 mL

■ VWR® UHP




2	1
0	0
4	0
2	0

20.42 μ L 1,000 μ L

Lock the volume

■ VWR® EHP

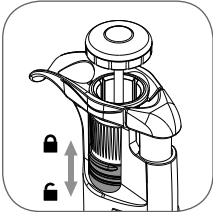


■ VWR® UHP





NOTE: VWR® Standard Single- and Multi-channel Pipettors do not feature a volume lock function.


Step 2

Insert the tip

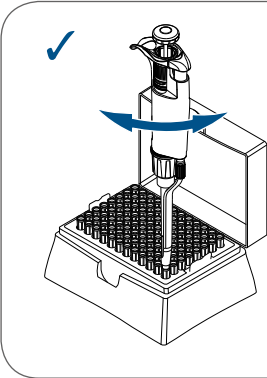

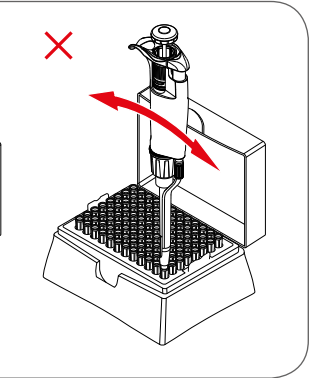


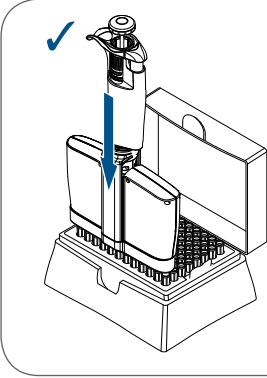

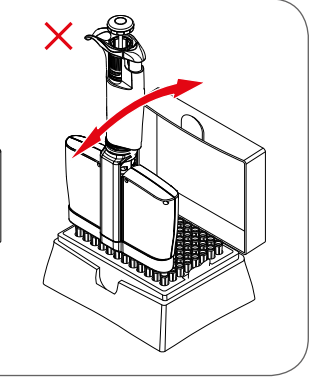


Ejector adjustment


VWR® Standard

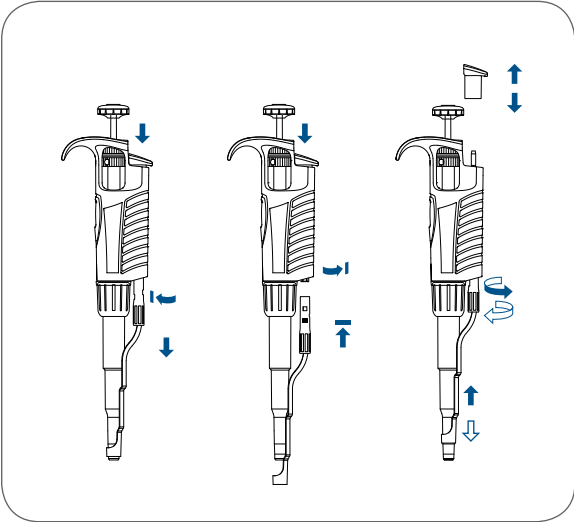

Pipettors 5,000 and 10,000 µL

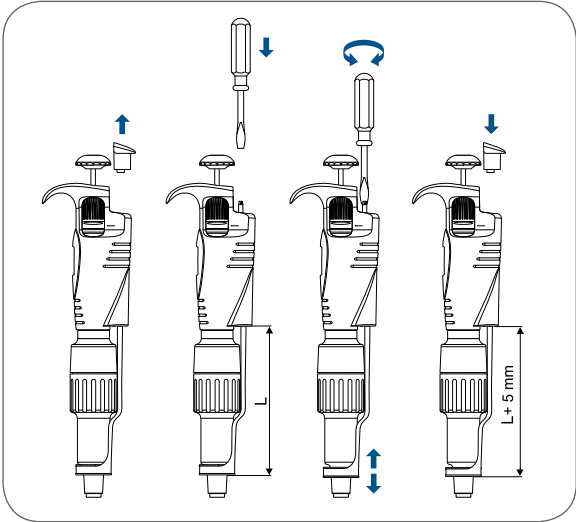



VWR® EHP

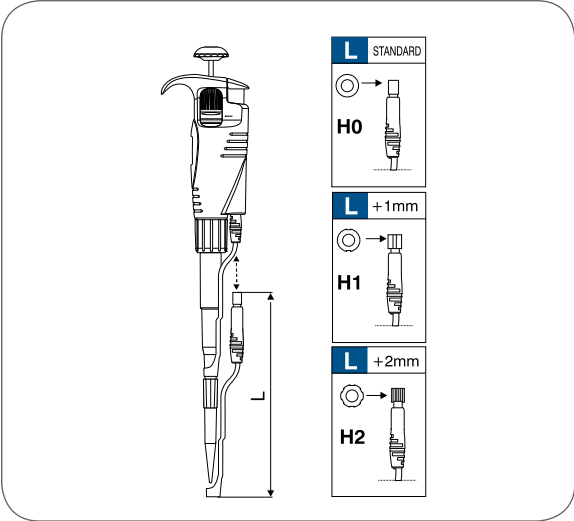


NOTE: Ejector adjustment is not necessary in VWR® UHP pipettors as they feature a long ejector stroke.

Step 3

Aspirate the liquid

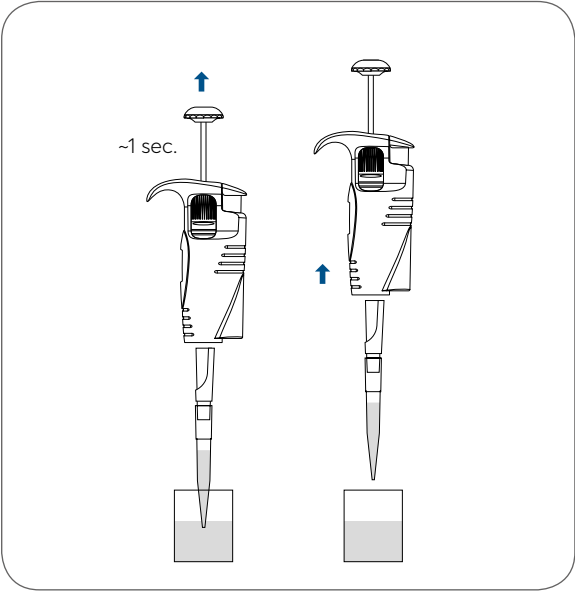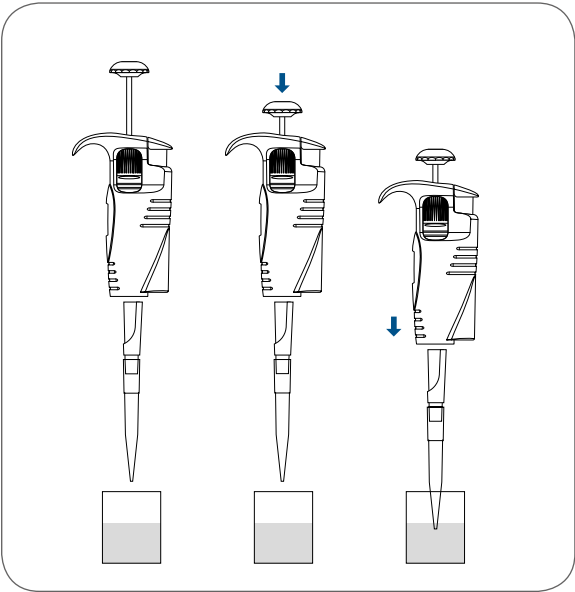


Recommended depth

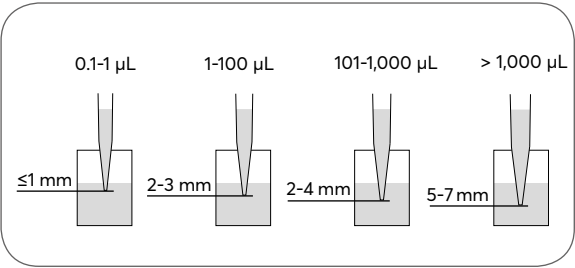


Step 4

Dispense the liquid

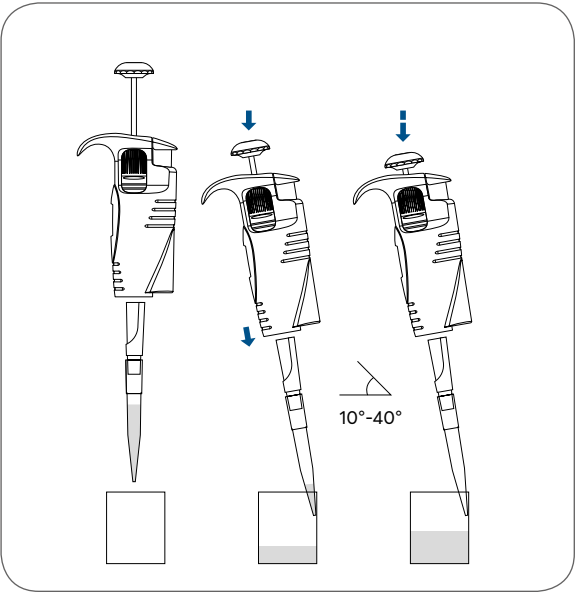

Step 5



Eject the tip

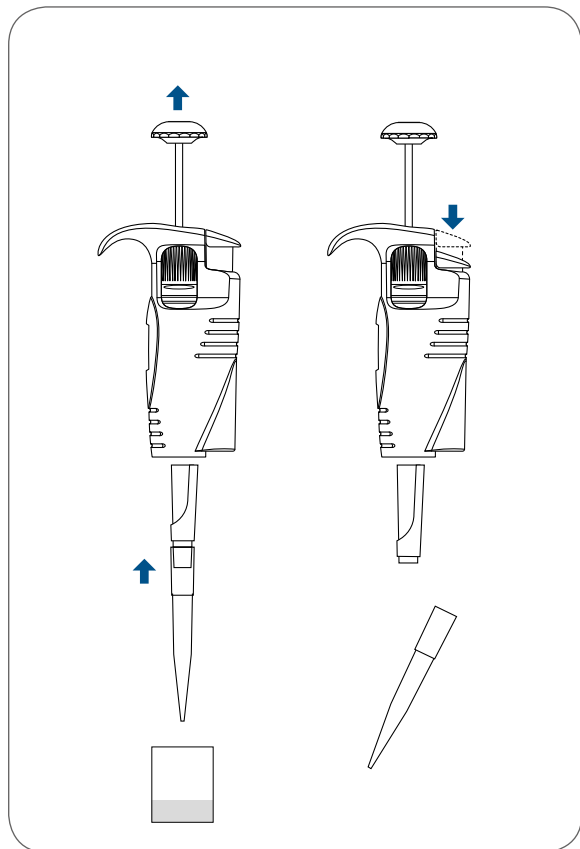

Step 6



Calibration

■ VWR® EHP ■ VWR® Standard

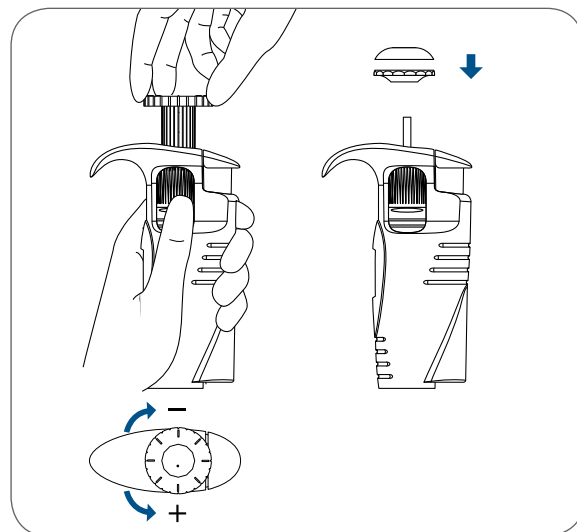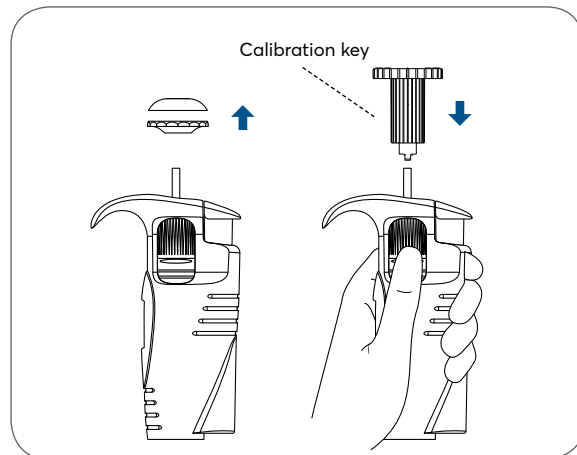

NOTE: Before calibration, in the VWR® EHP and VWR® UHP lines, always remember to lock the volume using the lock function and for the VWR® Standard hold the calibration knob.

■ VWR® UHP

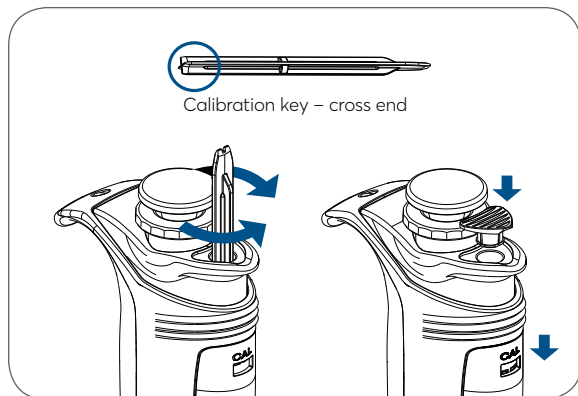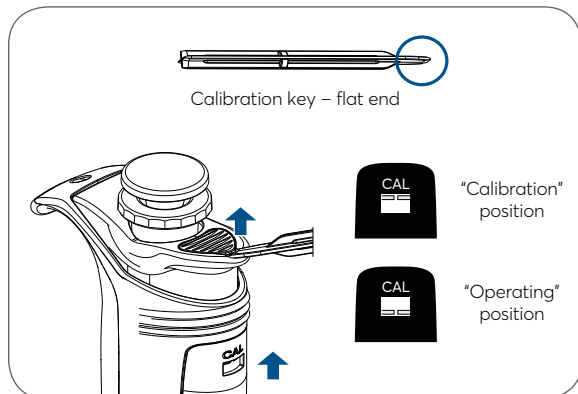

Warranty

VWR® Single- and Multi-channel Pipettors are covered by a three (3) year limited warranty. For more information on the warranty limitations refer to the full version of the User Manuals available at www.vwr.com

Safety Notes

- The pipettor is designed for the transfer of liquids only using the tip. Do not aspirate liquids without the tip attached. The aspirated liquid should not enter the pipettor, as it may cause damage.
- Single-use tips reduce the risk of contamination of samples.
- Keep the pipettor clean, avoiding the use of abrasive or corrosive cleaning agents (e.g., acetone).
- Keep the pipettor upright when there is liquid in the tip.
- Only using the pipettor in accordance with the manufacturer's instructions ensures the correct pipettor parameters are maintained.
- After replacing the plunger or the shaft, the pipettor should be calibrated.
- In the case of incorrect operation, the device should be cleaned in accordance with the Instructions for Use or transferred to a service point.
- Ambient operating temperature is +5°C to 45°C.
- Ambient storage conditions (in the original packaging during transport and short storage) is -25°C to 55°C.



- Follow general work safety regulations regarding hazards related to work in the laboratory.
- Take special care when pipetting aggressive substances.
- Use appropriate protective attire (e.g. clothing, goggles and gloves).
- Avoid pointing the pipettor at yourself or others during use.
- Only use parts and accessories recommended by the manufacturer.

Cleaning

- External parts may be cleaned with a swab moistened with isopropyl alcohol.
- Pipettors can be autoclaved at 121°C for 20 minutes. Prior to autoclaving, untighten the pipettor shaft, remove the filter from 5 and 10 mL pipettor shafts.
- The outer body of the pipettor is UV-resistant. The recommended distance from the radiation source to exposed element should not be less than 50 cm.
- Prolonged and very intense UV exposure can cause de-coloration of pipettor parts but does not affect its performance.

For additional product or technical information, visit www.vwr.com or contact your local sales office.

Austria

VWR International GmbH
Graumannsgasse 7
1150 Vienna
Tel.: +43 1 97 002 0
Email: info.at@vwr.com

Belgium

VWR International BV
Researchpark Haasrode 2020
Geldenaaksebaan 464
3001 Leuven
Tel.: +32 016 385 011
Email: vwr.be@vwr.com

China

VWR (Shanghai) Co., Ltd
Bld.No.1, No.3728 Jinke Rd,
Pudong New District
Shanghai, 201203- China
Tel.: +400 821 8006
Email: info_china@vwr.com

Czech Republic

VWR International s. r. o.
Veetee Business Park
Pražská 442
CZ - 281 67 Stránská Skalce
Tel.: +420 321 570 321
Email: info.cz@vwr.com

Denmark

VWR International A/S
Tobaksvej 21
2860 Søborg
Tel.: +45 43 86 87 88
Email: info.dk@vwr.com

Finland

VWR International Oy
Valimotie 17-19
00380 Helsinki
Tel.: +358 9 80 45 51
Email: info.fi@vwr.com

France

VWR International S.A.S.
Immeuble Estreo
1-3 rue d'Aurion
93114 Rosny-sous-Bois cedex
Tel.: 0 825 02 30 30* (national)
Tel.: +33 (0) 1 45 14 85 00
(international)
Email: info.fr@vwr.com
*0,18 € TTC/min + prix appel

Germany

VWR International GmbH
Hilpertstrasse 20a
D - 64295 Darmstadt
Freecall: 0800 702 00 07
Tel.: +49 (0) 6151 3972 0
(international)
Email: info.de@vwr.com

Hungary

VWR International Kft.
Simon László u. 4.
4034 Debrecen
Tel.: +36 (52) 521-130
Email: info.hu@vwr.com

India

VWR Lab Products Private
Limited
No.139, BDA Industrial Suburb,
6th Main, Tumkur Road,
Peenya Post,
Bangalore, India - 560058
Tel.: +91-80-28078400
Email: vwr_india@vwr.com

Ireland

Northern Ireland

VWR International Ltd /
VWR International (Northern
Ireland) Ltd
Orion Business Campus
Northwest Business Park
Ballycoolin
Dublin 15
Tel.: +353 01 88 22 222
Email: sales.ie@vwr.com

Italy

VWR International S.r.l.
Via San Giusto 85
20153 Milano (MI)
Tel.: +39 02-3320311/02-487791
Email: info.it@vwr.com

Korea

VWR International -
17, Daehak 4-ro, Yeongtong-gu,
Suwon-si, Gyeonggi-do
Tel.: +82 31 645 7256
Email: saleskorea@avantorsciences.com

Middle East & Africa

VWR International FZ-LLC
DSP Laboratory Complex
125, Floor 01
Dubai, United Arab Emirates
Tel.: +971 4 5573271
Email: info.mea@vwr.com

The Netherlands

VWR International B.V.
Postbus 8198
1005 AD Amsterdam
Tel.: +31 020 4808 400
Email: info.nl@vwr.com

Norway

VWR International AS
Brynsalleen 4
0667 Oslo
Tel.: +47 22 90 00 00
Email: info.no@vwr.com

Poland

VWR International Sp. z o.o.
Limbowa 5
80-175 Gdansk
Tel.: +48 058 32 38 210
Email: info.pl@vwr.com

Portugal

VWR International - Material
de Laboratório, Lda
Centro Empresarial de
Alfragide
Rua da Indústria, nº6
2610-088 Amadora
Tel.: +351 21 3600 770
Email: info.pt@vwr.com

Singapore

VWR Singapore Pte Ltd
18 Gul Drive
Singapore 629468
Tel.: +65 6505 0760
Email: sales.sg@vwr.com

Spain

VWR International Eurolab S.L.
C/ Tecnología 5-17
A-7 Llinars Park
08450 - Llinars del Vallès
Barcelona
Tel.: +34 902 222 897
Email: info.es@vwr.com

Sweden

VWR International AB
Fagerstagatan 18a
163 94 Stockholm
Tel.: +46 08 621 34 00
Email: info.se@vwr.com

Switzerland

VWR International GmbH
Lerzenstrasse 16/18
8953 Dietikon
Tel.: +41 044 745 13 13
Email: info.ch@vwr.com

UK

VWR International Ltd
Customer Service Centre
Hunter Boulevard
Magna Park
Lutterworth
Leicestershire
LE17 4XN
Tel.: +44 (0) 800 22 33 44
Email: uksales@vwr.com

United States

VWR International LLC
100 Matsonford Rd
Radnor, PA 19087
+1 800-932-5000
www.vwr.com

Manufactured for VWR International, LLC
100 Matsonford Rd, Radnor, PA 19087

Version 1

Made in Poland